PERISTA® Pump

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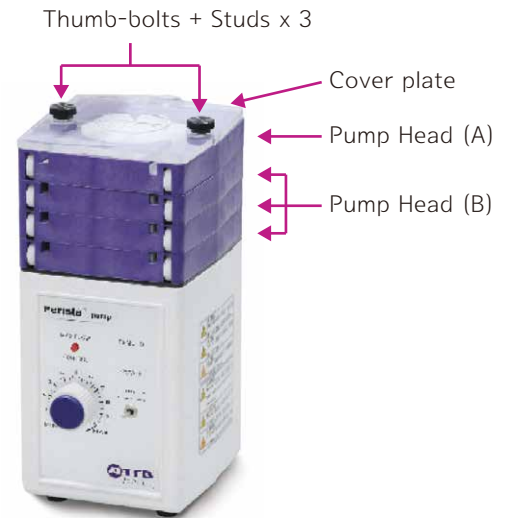
Pump Head (A) : 최상단 pump head

Ordering Information

Cat. No.	Product Name
1292231	Pump Head (A) * Rollers included, Cover plate Not included
1292232	Pump Head (B) * Rollers included, Cover plate Not included
1292233	Set of rollers (A)
1292234	Set of rollers (B)
1292235	Cover plate
1292236	Addition kit [Pump head(B) + 2 Studs + silicon tube(1m each)]
1292133	Thumb-bolts 2/pk
1292213	Studs 2/pk

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O.D. : outside diameter (외경)
I.D. : inside diameter (내경)

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Silicon Tube Chemical Proof Chart

◎ Excellent, almost completely unaffected
○ Slightly affected but presents no problem for usage

△ Infiltrated to a certain degree, not recommended
× Extreme damage not usable

Sodium nitrite	◎	Ricinus	◎	Phenol	○	Potassium hydroxide	◎	Hydrogen peroxide solution	○
Asphalt	×	Sodium hydroxide	○	Butadiene	×	Chloroform	×	Ammonium hydrogen carbonate	◎
Acetaldehyde	×	Paraldehyde	×	Butanol	△	Chlorobenzene	×	Aromatic carbon hydride	×
Acetone	○	Formic acid	◎	Butane	×	Chloromethane	×	Chromic acid anhydride	◎
Aniline	×	Styrene	×	Fulfural	×	Kerosene	×	Ammonium hydroxide	◎
Linseed oil	◎	Stearic acid	△	2-propanol	○	Magnesium hydroxide	◎	0~10% sulfuric acid	◎
2-aminoethanol	○	Fat acid	○	Benzene	×	Trichloroacetic acid	△	10~50% sulfuric acid	△
Sulfurous acid	△	Tannic acid	◎	Pentanol	○	Acetic acid	○	Methyl isobutyl ketone	×
Sodium sulfite	◎	Diesel oil	○	Pentane	×	Acetic ether	×	Acetic acid anhydride	○
Ammonia	○	Molasses	◎	Benzene	×	Sodium acetate	○	Phtanol acidanhydride	○
Ammonia water	○	Triethanolamin	◎	Maleic acid	◎	Butyl acetate	×	Methyl ethyl ketone	×
Isobutane	×	Triethylamine	◎	Palmitic acid	○	Propyl acetate	×	Monochloroacetic acid	○
Ethanol	○	Triol	×	Silicon oil	△	Pentyl acetate	×	50% sulfuric acid	×
Ethylene glycol	○	Trichloroethane	×	Pitch	×	Methyl acetate	×	2-methyl propene	×
Hydrochloric acid	×	Coal tar	△	Methanol	◎	Calcium hypochlorite	◎	Chlorine dielectric oil	○
Xylene	×	Naphtha	×	Nitric acid	×	Sodium hypochlorite	△	Chlorine (moistened)	◎
Fish oil	△	Sulfur dioxide	△	Creosote	×	Diamylamine	◎	Dimethyl phthalate	△
Octanol	◎	Nitrobenzene	×	Cotton oil	◎	1,2-dichloroethane	×	Paraform aldehyde	×
Oleic acid	○	Lactic acid	◎	Melamine resin	○	Dichloropentane	×	Kalium phosphate	△
Glycerol	◎	Hydroquinone	○	Melt sulfur	△	Dipentane	×	10% sodium hydroxide	◎

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Providing research solutions for
Biochemistry / Molecular Biology /
Genetic engineering development

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